


A-0214 Vacu-vials® Snapper

(An alternative to the sample cup for use with K-5543 Peroxide and K-7433 Ozone Vacu-vials ampoules)

1. Place the Vacu-vial ampoule, tip first, into the snapper using care not to prematurely snap the tip of the ampoule.
2. Place the snapper with ampoule into the sample to be tested.

Note: There must be enough water in the sample container to cover at least the bottom half of the snapper.

3. Apply gentle downward pressure to the bottom of the ampoule to snap the tip. The ampoule will fill leaving a bubble for mixing.
4. Complete the specific test procedure provided in the Vacu-vials kit being used.

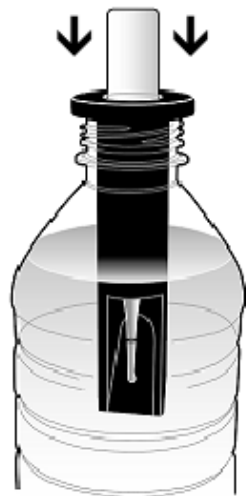

www.chemetrics.com
4295 Catlett Road, Midland, VA 22728 U.S.A.
Phone: (800) 356-3072; Fax: (540) 788-4856
E-Mail: orders@chemetrics.com
Feb. 2022, Rev. 2