A-0214 Vacu-vials® Snapper

(An alternative to the sample cup for use with K-5543 Peroxide and K-7433 Ozone Vacu-vials ampoules)

- 1. Place the Vacu-vial ampoule, tip first, into the snapper using care not to prematurely snap the tip of the ampoule.
- Place the snapper with ampoule into the sample to be tested.

Note: There must be enough water in the sample container to cover at least the bottom half of the snapper.

- Apply gentle downward pressure to the bottom of the ampoule to snap the tip. The ampoule will fill leaving a bubble for mixing.
- 4. Complete the specific test procedure provided in the Vacu-vials kit being used.



www.chemetrics.com 4295 Catlett Road, Midland, VA 22728 U.S.A. Phone: (800) 356-3072; Fax: (540) 788-4856 E-Mail: orders@chemetrics.com Apr. 2023, Rev. 3